

PBST Buffer (20X)

Catalog number: ACC 00501

Lot number	Item
_____	50 ml - makes 1 liter

Store PBST buffer (20X) concentrate at room temperature for up to one year.

Preparing Buffer from concentrate

PBST buffer is used in all Agdia ELISA wash steps. Additionally it can be used as a sample extraction buffer in some tests and also as a base for many other buffer formulations.

Dilute 1 volume of 20X PBST buffer concentrate with 19 volumes of distilled water before use.

Example: To prepare 1 liter of working 1X PBST buffer

Buffer concentrate	50 ml
Distilled water	950 ml

Measure the distilled water into a container and add the buffer concentrate while stirring.

Storage

Agdia recommends preparing only as much 1X PBST buffer as is needed for one day.