



**BioFX® Alkaline Phosphatase/Horseradish
Peroxidase Block**
Product insert

Product number:

APHP-0100-01 (100 mL)

Intended use:

BioFX® Alkaline Phosphatase/Horseradish Peroxidase Block is supplied as a ready-to-use blocking reagent for immunohistochemistry samples. The solution blocks both alkaline phosphatase and peroxidase endogenous enzymatic activity.

Product stability and storage:

Product stability	3 years from the date of manufacture.
Storage	Store at 2-8°C. Care should be taken to minimize the risk of contamination.

Recommendations for use:

- 1) Samples should be blocked for a minimum of 10 minutes at room temperature (25°C).
 - Additional blocking time may be required depending on tissue thickness and sample penetrability.
- 2) Samples should be rinsed with copious amounts of distilled water to remove excess blocking solution.

Additional considerations:

For technical assistance, email ivdtechsupport@surmodics.com

Related Products:

BioFX® DAB Immunohistochemistry Substrate (DABI)

BioFX® Fluorescence Preserving Media (FLPM)

All products listed are for research use or for further manufacturing into *in vitro* diagnostic reagents. The products are not intended for use in humans or animals. Sales are without any seller's warranty or representation, expressed or implied, by usage or otherwise; no claims beyond replacement of product not meeting specifications or refund of purchase price shall be allowed. All claims must be made within 30 days following date of delivery.