

user instructions



711 HULMAN STREET
TERRE HAUTE, IN 47802
812-235-6167
FAX 812-234-6975

HOTSPOT Heat Lamp

108A HS250-3



The HOTSPOT Heat Lamp is shipped assembled. The 250W infrared lamp is screwed in its socket in the housing and is protected by cushioning material. Remove and discard this material before use.

To facilitate packaging and shipping, the one-half-inch diameter mounting rod is folded back against the housing. For use, bring it into the desired position by simply exerting enough force to overcome the resistance of the spring-loaded mounting joint. When you clamp the rod on an apparatus frame or ring stand, you can easily aim it in the desired direction. The high quality ball-and-socket joint is made to hold the lamp firmly in position and give years of service without requiring adjustment.

To operate the device, simply plug the cord into a 120V outlet and turn on the switch. On variable control models, the switch is activated by pushing down on the Control Knob. After the lamp is turned on, rotating the Knob clockwise will increase the voltage to give the desired intensity of heat. Caution: The housing of a heat lamp gets hot after a period of use. Use care when redirecting the position of the lamp when it is ON, or has just been turned OFF.

LIMITED WARRANTY

Glas-Col warrants products of its manufacture to be free from defects in materials and workmanship for 1 year from date of purchase. We will repair or replace without charge any product found to be defective {upon examination at the factory} within this one year period, providing that the apparatus has not been overheated, abused, misused, overloaded or subjected to spillage. Glas-Col is not responsible for damages to apparatus due to improper installation or through attempts to operate the apparatus beyond its rated capacity. Normally expendable parts are not covered by this warranty.

LIMITATION OF WARRANTIES

APART FROM SUCH WRITTEN STATEMENT OF WARRANTY, THERE ARE NO WARRANTIES, EXPRESSED, IMPLIED OR STATUTORY, WHICH EXTEND BEYOND THE DESCRIPTION OF THE PRODUCTS ON THE FACE HEREOF.

Glas-Col products are intended only for legal and legitimate purposes in commercial laboratory settings.

Glas-Col reserves the right to make product refinements without prior notice.

Glas-Col,LLC

711 Hulman Street., P.O. Box 2128, Terre Haute, IN 47802-0128

Phone. (812) 235-6167, Fax. (812) 234-6975

e-mail: pinnacle@glascol.cpm

Web: www.glascol.com