



## **Instructions for Washing Gerber Equipment**

**The following glass cleaning procedures are recommended:**

Laboratory glassware with graduations and imprints can be cleaned manually or mechanically. However, care must be taken to ensure that the rinsing bath temperature does not exceed 70 °C when using alkaline detergents.

In addition, the time spent in the rinsing bath must be kept to a minimum. Alkaline media in combination with temperatures above 50 °C and longer residence times can lead to volume changes due to glass removal and to the destruction of the prints.

Our recommendation for cleaning is water above 90 °C without chemical additives (hot water cleaning).