

# user instructions



**Glas-Col®**

711 HULMAN STREET  
PO BOX 2128  
TERRE HAUTE, IN 47802  
812-235-6167  
FAX 812-234-6975

## **LABORATORY STAND MODEL NO. L70**

The L70 laboratory stand is designed to help support Glas-Col stirrer motors and homogenizing equipment. The L70 can also be used for general laboratory use. The L70 provides a sturdy horseshoe shaped base with four feet. A 23" long, 3/4" diameter vertical support rod attaches to the base by means of a thumb screw.

### **PARTS INCLUDED:**

- 1 - Base with Rubber Bumper Feet (R68 109)
- 1 - 3/4" Dia. Rod
- 1 - 5/16-18 x 1" Thumb Screw

### **ASSEMBLY:**

Place the 3/4" diameter support rod down through the hole in the base. Make sure flat on support rod is aligned with the tapped hole in the base and tighten thumb screw.

### **MAINTENANCE:**

The lab stand should be protected from spills, damage and corrosive atmospheres so far as possible. Any spills should be cleaned quickly and any damage should be repaired immediately. In time it may be necessary to replace some parts of the lab stand. A list of common replacement parts are listed below:

**Rubber Bumper Feet.....099B R68 109**  
**5/16-18 x 1" Thumb Screw.....099B R88 208**

### **Limited Warranty**

Glas-Col warrants products of its manufacture to be free from defects in material and workmanship for one year and agrees to repair or replace without charge any products found defective upon examination at the factory. With proper care and operation, Glas-Col products will give long and efficient service. Chemical spillage, overheating, overloading, and general misuse will greatly reduce the service life. Glas-Col is not responsible for damage to apparatus due to improper installation or through attempts to operate the apparatus beyond its rated capacity, intentional or otherwise. Normally expendable parts are not covered by this warranty.

### **LIMITATION OF WARRANTIES**

**APART FROM SUCH WRITTEN STATEMENT OF WARRANTY, THERE ARE NO WARRANTIES, EXPRESSED, IMPLIED OR STATUTORY WHICH EXTEND BEYOND THE DESCRIPTION OF THE PRODUCTS ON THE FACE HEREOF.**

Glas-Col products are intended only for legal and legitimate purposes in commercial laboratory and industrial settings.

Glas-Col reserves the right to make product refinements without prior notice.

#### **Glas-Col, LLC**

711 Hulman Street, P.O. Box 2128, Terre Haute, IN 47802-0128

Phone. (812) 235-6167, Fax. (812) 234-6975

E-mail: [pinnacle@glascol.com](mailto:pinnacle@glascol.com)

Web: [www.glascol.com](http://www.glascol.com)